

OCEAN OPTICS UV/VIS SPECTROMETERS INSTRUCTIONS FOR START-UP/SHUT

DOWNÍ Dell Computers

Updated November 14, 2017

Celeste Morin, x3756, Druckenmiller 243

- 1) Plug the power cable into the wall. The spectrometer fan may turn on at this point. If not, it will turn on in Step #2.
- 2) Plug the USB cable into the spectrometer and USB port in the front of the computer. Fan should be on at this point, if not see your lab instructor.
- 3) Open the Spectrasuite software (rainbow-colored icon). Communication has been established if you see a picture of the UV-VIS spectrometer in the main window of the software.
- 4) When finished using the spectrometer, close the Spectrasuite software, unplug the USB cable from the computer, and unplug the power cable from the wall.